

CERTOCLAV VACUUM PRO

THE LABORATORY STERILIZER WITH A REVOLUTIONARY
USER INTERFACE



**AUTOMATIC
DOCUMENTATION
IN THE CLOUD**


**ON MACHINE
USER SUPPORT**

HIGHLY DEVELOPED FEATURES FOR COMFORTABLE USE




Mobility

- Mobile autoclave display can be placed at your favorite position.
- Included mountable industrial stand ensures save stand.

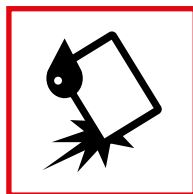


Durability

- Mobile display with certified industrial grade.
- IP 64 protection and shock resistant.

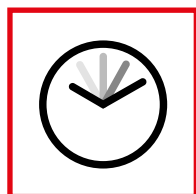

High Performance

- Sterilisation of solid instruments within 20 minutes (Time including fractionated pre-vacuum and drying).

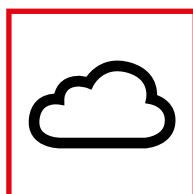

Innovation

- Connection of multiple autoclaves to local network.
- CertoCloud via Ethernet cable or wireless LAN.



ACCESSORY

Tablets

Art.-No.:

CV-8-T-S-Pro
CV-12-T-S-Pro
CV-18-T-S-Pro
CV-22-T-S-Pro



Extra set of tablets for all models of CertoClav Class B sterilizers.

Water Distiller

Art.-No.:

CV-WD-3L
CV-WD-5L

With a volume of 3 or 5 liters.



Sealing machine

Art.-No.:

CV-SM-M
CV-SM-L


2 sizes, incl. different types of packing strap.



CertoCloud

EASY DATA ACCESS AND STORAGE. ONLINE DEVICE MANAGEMENT.

- Any kind of documentation according to the customers needs.
- Logging via USB-Stick, CF-Card, direct printer output or fully automatic to the CertoCloud.
- Accessible via www.certocloud.com or as offline application.

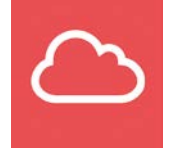

Secure Connection

- SSL encryption
- Secure server
- 100% data privacy




Multi Device Management

- Combine all devices in one account
- Create programs
- Disseminate programs to all devices



Long Term Data Storage

- All data stored for 10 years to fulfil documentation requirements
- Secure backup

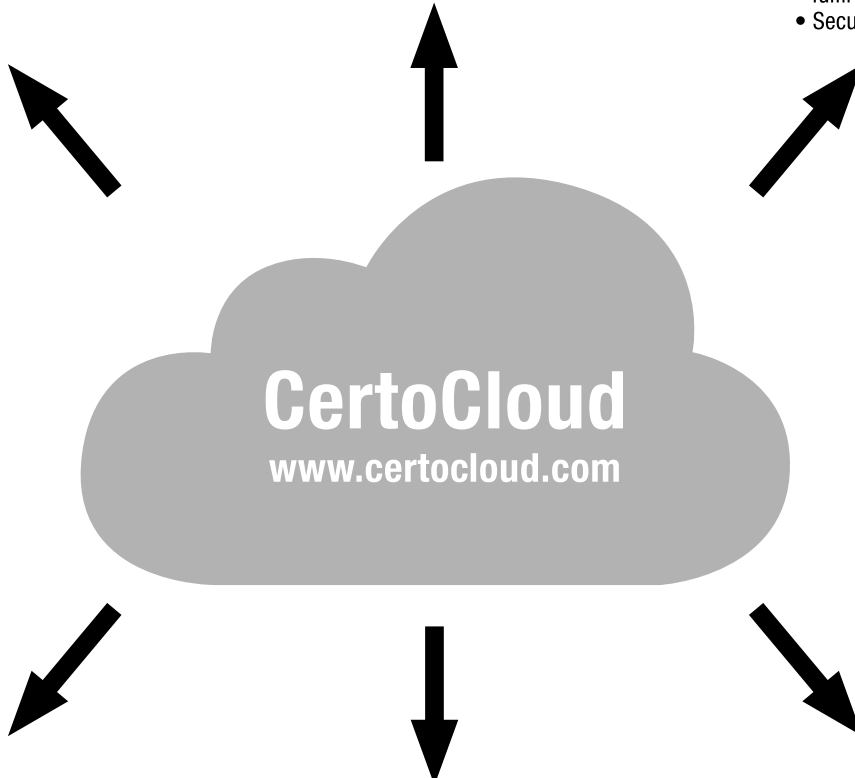

Customization

- Custom reports
- Direct printing
- Freely definable data formats



Email Notifications

- Alarms
- Cycle complete
- Process data



User Management


- Individual accounts
- Accounts linked to devices

THE USER INTERFACE WITH ON-SCREEN TRAINING.



Step-by-Step video guides

- All usage steps explained
- No more need to train staff
- Videos can be skipped
- Ensures simple usage

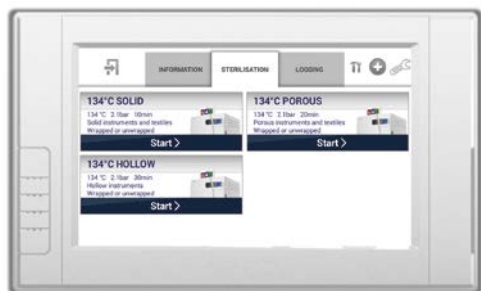

Program Selection


- Detailed description
- Custom program

PROGRAMS	Vacuum Times	Sterilizing Time	Sterilizing Temperature	Sterilizing Pressure	Vacuum Persist Time	Dry Time
121° C SOLID	1	20 min	121° C	110 kPa		3 min
121° C POROUS	3	20 min	121° C	110 kPa		7 min
121° C HOLLOW	3	20 min	121° C	110 kPa		10 min
134° C SOLID	1	4 min	134° C	210 kPa		3 min
134° C POROUS	3	4 min	134° C	210 kPa		7 min
134° C HOLLOW	3	4 min	134° C	210 kPa		10 min
USER DEFINED	3	5 min	134° C	210 kPa		10 min
B-D TEST	3	3,5 min	134° C	210 kPa		7 min
VACUUM TEST				-80 kPa	15 min	
CLEAN PROCESS	3	5 min	105° C	20 kPa		10 min
PRIONEN	3	19 min	135° C	210 kPa		10 min




Protocol history

- Protocols stored locally
- Graphically displayed
- Errors explained detailed

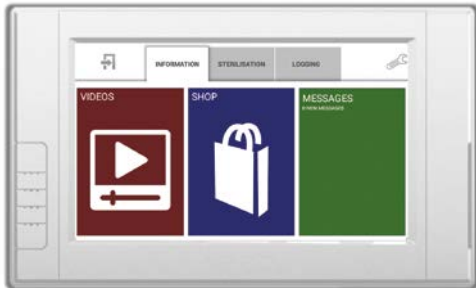

Service

- Order spare parts On-Screen
- Message center
- CertoClav Academy videos

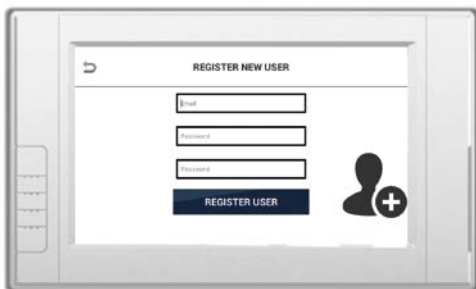

Advanced Multi-User Management

- Secure user management system that enable numerous members of personnel.
- Personal email and password, can also be used for CertoCloud account.
- User defined settings and accurate association of performed cycles with operators.



TECHNICAL DETAILS

	Vacuum Pro 8 L	Vacuum Pro 12 L	Vacuum Pro 18 L	Vacuum Pro 22 L
Art. Nr.	CV-8-Vac-Pro	CV-12-Vac-Pro	CV-18-Vac-Pro	CV-22-Vac-Pro
Description	Vacuum autoclave, Class B, GLP ready, Stand-alone			
Compliance	CE			
Rated Voltage, Output	220V, 50 Hz, 2400W			
Temperature Range	121° C - 134° C			
Max working pressure	0,23 MPA			
Chamber material	Stainless steel 304			
Interface	7 inch touch screen			
Safety	Electrical locker			
Equipment	Optional label printer, USB-Stick, Ethernet, Wireless LAN, CF-Card			
Product size	53,5 x 36,5 x 32 cm	53,5 x 36,5 x 32 cm	60 x 50 x 43 cm	60 x 50 x 43 cm
Diameter	17 cm	20 cm	25 cm	25 cm
Depth	32 cm	36 cm	35 cm	45 cm
Volume	8 l	12 l	18 l	22 l
Weight	52 kg	52 kg	65 kg	65 kg
Languages	English, German			
CertoCloud Licence	Unlimited			
Warranty	2 years			

CERTOCLAV

Certoclav Sterilizer GmbH
 Georg Grinninger-Str. 37
 A-4050 Traun
 Phone: +43 (0)7229 689
 www.certoclav.com

Orders, technical customer service, custom-manufacture, dealer verification:

Christina Bokesch, Phone: +43 (0)7229 689 26, Fax: +43 (0)7229 74 141
 Email: bokesch@certoclav.com

Product Manager: Klaus David, Phone: +43 (0)7229 689 28, Email: david@certoclav.com

Operating instructions, technical information and operation videos: www.certoclav.com.